


# Powdermax®

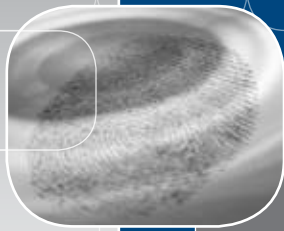

Esco Powdermax 1 Powder Weighing Balance Enclosure  
Model PW1-3A\_

## Powder Weighing Balance Enclosure

Laboratory operations involving weighing, mixing and dispensing hazardous or nuisance powders can now be performed safely and economically with the Esco Powdermax 1 Powder Weighing Balance Enclosure. The enclosure's unique design maintains balance stability to 4 decimal places without compromising containment.

Designed specifically for powder containment applications, the Powdermax Enclosure employs carefully directed inward airflow at the face of the enclosure to protect the operator from exposure to hazardous powders or potent pharmaceutical compounds. The aerodynamically designed enclosure reduces airflow turbulence for maximum containment while minimizing the effect that air drafts may have on sensitive weighing balances. The enclosure design exceeds the containment requirements of ANSI/ASHRAE 110 as proven by independent type testing. All exhaust air is HEPA-filtered, removing virtually all particulate matter, before recirculation to the laboratory.

### Designed for Enhanced Usability and Efficiency

The Powdermax 1 is designed specifically to provide the operator with a high level of usability, comfort and visibility.


- A 13-degree sloped front allows easy access to the work zone, eliminating operator fatigue and increasing productivity.
- Transparent 6 mm (0.2") frameless acrylic front and sides enhance visibility and operator comfort during longer periods of operation. The self-supporting pre-tensioned hinges on the front window provide easy access during loading and start-up.
- Ergonomically designed apertures for hands simultaneously provide maximum movement within the work zone while protecting the operator. The curved front edge minimizes airflow turbulence and improves user comfort.

- Electronic ballast for the fluorescent lighting provides zero-flicker with increased energy efficiency, reliability and service life with a lower heat output.
- Two electrical pass throughs on the back wall of the unit provide convenient access to power sources for equipment.
- The spacious interior dimensions will easily accommodate large analytical balances.

### Enhanced Filtration System

The enclosure's unique filtration system ensures the highest level of protection for the operator.

- The Powdermax 1 provides enhanced safety to laboratory personnel by drawing air across the front opening in the hood at a controlled rate, preventing the escape of powders and particulate matters into the general laboratory environment.
- A unique filter mounting system eliminates bypass leaks inherent in conventional clamping systems.
- The Powdermax 1 is equipped with a generously sized HEPA filter system, rated at >99.99% efficiency for 0.3 micron sized particles.
- The aerodynamically designed system constructed of 1.2mm epoxy powder-coated electrogalvanized steel delivers maximum containment by ensuring airflow uniformity throughout the work chamber of the cabinet.
- Since HEPA-filtered air is recirculated to the laboratory, the enclosure does not have to be connected to an extraction/exhaust system, resulting in low operating costs and mobility.

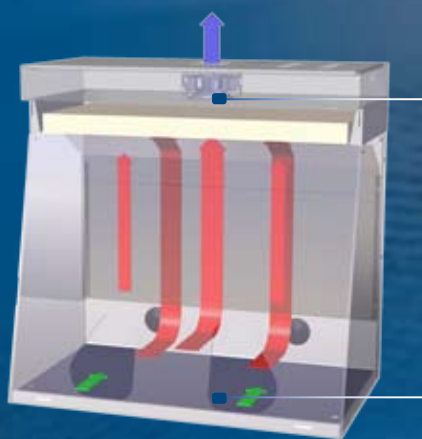

## Powdermax 1 Powder Weighing Balance Enclosure Airflow

Blower  
Supply HEPA Filter

- HEPA-filtered air
- Unfiltered / potentially contaminated air
- Inflow air

- Room air is drawn in across the front of the cabinet with an average velocity of 0.5 m/s (100 fpm).
- Air is drawn up through the cabinet's work zone and forced through the HEPA filter (>99.99% typical efficiency for 0.3 micron sized particles).
- The state-of-the-art baffle system ensures airflow uniformity throughout the cabinet's main chamber.
- The HEPA filtered air then returns to the laboratory stripped of all airborne contaminants or is vented through the optional exhaust collar to exhaust ducting for enhanced safety.

### User-Friendly Control System

The user-friendly Esco Sentinel™ micro-processor based control system is fully configurable according to operator's requirements and comes equipped with a number of enhanced features to promote cabinet usability.

- Accurate true airflow velocity sensing technology measures all critical cabinet airflow parameters allowing superior monitoring. Temperature compensated sensors ensure increased accuracy.
- Password-protected administration can be set to restrict access to the main menu.
- Solid state variable speed controllers offer superior control over conventional "step" controllers. The built in RFI and noise filters eliminate interference with adjacent instrumentation.
- A bright, easy-to-read, LCD display provides continuous monitoring of cabinet airflow.
- Audible and visual alarms ensure operator and environmental protection by alerting the user in the event of low airflow or unsafe sash positions.
- When the cabinet is switched on, its automatic warm-up cycle is activated, purging all contaminants from the work zone prior to use.

- An automatic post-purge cycle can be configured on shutdown, purging all residual contaminants before deactivation.

### Highest Quality Enclosure Construction

All Esco cabinets are constructed to the highest quality using the finest materials.

- The enclosure's head unit is constructed of electrogalvanized steel with abrasion-resistant powder-coated finish.
- External surfaces are coated with Esco Isocide™ antimicrobial coating to protect against surface contamination and inhibit bacterial growth. Isocide eliminates 99.9% of surface bacteria within 24 hours of exposure.
- The enclosure is supplied with a chemical and abrasion resistant phenolic resin removable work surface. The solid work surface minimizes transmission of fan vibration to the user's weighing balance.

### Blower Efficiency

The Powdermax 1 incorporates an energy efficient external rotor motor design which reduces operating costs and has extremely low noise and vibration levels. The cabinet blower system creates an inflow of air from

the ambient environment into the cabinet with an average velocity of 0.4 m/s (80 fpm).

### Designed and Built to Exceed Safety Criteria

All components used in Esco products meet or exceed all the applicable safety requirements and are UL listed/recognized. The Powdermax 1 is compliant with International Standards. The cabinet is designed to meet the general safety requirements of EN 61010-1, EN 61326 Class B.

### Options and Accessories

The Powdermax 1 is available with a number of options and accessories to meet your needs. These include:

- Mobile cart with foldable trays
- Height-adjustable lab chair
- Ergonomic foot rest

Further customization specifications are available upon request.

### Warranty

The Powdermax 1 is warranted for 3 years excluding consumable parts and accessories. Contact your local sales representative for specific warranty details.

	Containment	Electrical Safety
Standards Compliance	ANSI/ASHRAE 110-1995, USA	UL-61010-1, USA CAN/CSA-C22-2 No.61010-1 EN-61010-1, Europe IEC61010-1, Worldwide

## Powdermax®

Specialty Cabinets • Powder Weighing Balance Enclosures


## General Specifications, Powder Weighing Balance Enclosure, Model\_ PW1-3A

*\*Note to Customer: Insert electrical voltage number into last number digits \_ when ordering.*

Nominal Size	0.9 meters (3 ft)	
External Dimensions (W x D x H)	840 x 700 x 1125 mm (33" x 27.6" x 44.3")	
Internal Work Area, Dimensions (W x D x H)	820 x 650 x 848 mm (32.3" x 25.6" x 33.5")	
Usable Work Zone	0.52 m <sup>2</sup> (5.6 sq.ft)	
Air Volume (at Initial Velocity)	234 m <sup>3</sup> /h (138 cfm)	
Inflow Velocity	Initial setpoint: average of 0.5 m/s (100 fpm) measured in plane of work aperture	
HEPA Filter Efficiency	>99.99% at 0.3 microns	
Sound Emission per NSF / ANSI 49	<55 dBA at initial blower speed setting measured at typical operator work position	
Fluorescent Lamp Intensity at Zero Ambient	>560 Lux (> 52 foot-candles)	
Cabinet Construction	Main Body	1.2 mm / 0.06" / 16 gauge electro-galvanised steel with white oven-baked epoxy powder-coated finish.
	Work Zone	6 mm (0.2") Acrylic
	Side Walls	6 mm (0.2") Acrylic
Power Rating	225W (230V, AC units) 270W (115V, AC units)	
Net Weight **	66 kg (145.5 lbs)	
Shipping Weight **	96.4 kg (212.5 lbs)	
Shipping Dimensions, Maximum (W x D x H)**	1000 x 1100 x 1180 mm (39.8" x 43.3" x 46.5")	
Shipping Volume, Maximum**	1.3 m <sup>3</sup> (45.9 cu.ft.)	
Electrical*	<b>Model</b>	<b>Voltage</b>
	PW1 - 3A1	220-240V, AC, 50Hz, 1Ø
	PW1 - 3A2	110-120V, AC, 60Hz, 1Ø
	PW1 - 3A3	220-240V, AC, 60Hz, 1Ø

\* Additional voltages may be available; contact Esco for ordering information.  
 \*\* Cabinet only; excludes optional stand.

### Powdermax 1 Powder Weighing Balance Enclosure Technical Specifications

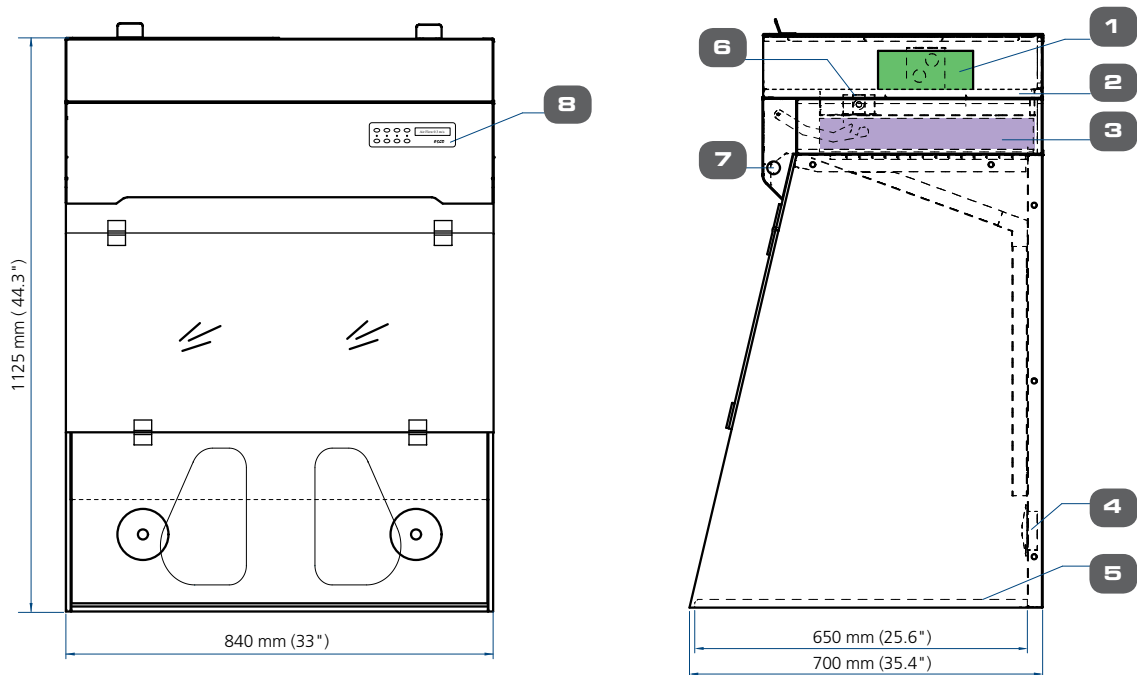

- |                     |                               |                      |                                    |
|---------------------|-------------------------------|----------------------|------------------------------------|
| 1. Blower           | 3. HEPA Filter                | 5. Phenolic Work Top | 7. Fluorescent Light               |
| 2. Electrical Panel | 4. Service Fixture Hole Cover | 6. Sensor Box        | 8. Sentinel Microprocessor Control |



**Esco Containment, Clean Air and Laboratory Equipment Products**

- Biological Safety Cabinets, Class II, III
- Fume Hoods, Conventional, High Performance, Ductless Carbon Filtered
- Laminar Flow Cabinets, Horizontal, Vertical, PCR
- Animal Containment Workstations
- Hospital Pharmacy Isolators, Cytotoxic Safety Cabinets
- Specialty Workstations: *In-Vitro* Fertilization, Powder Weighing
- PCR Thermal Cyclers, Conventional, Real-Time
- Cleanroom Fan Filter Units, Modular Rooms, Air Showers, Pass Thrus

Since 1978, Esco has emerged as a leader in the development of controlled environment, laboratory and cleanroom equipment solutions. Products sold in more than 100 countries include biological safety cabinets, fume hoods, ductless fume hoods, laminar flow clean benches, animal containment workstations, cytotoxic cabinets, hospital pharmacy isolators, and PCR cabinets and instrumentation. With the most extensive product line in the industry, Esco has passed more tests, in more languages, for more certifications, throughout more countries than any biosafety cabinet manufacturer in the world. Esco remains dedicated to delivering innovative solutions for the clinical, life science, research and industrial laboratory community. [www.escoglobal.com](http://www.escoglobal.com).

NSF / ANSI 49 Biological Safety Cabinets • Animal Containment Workstations • Fume Hoods • Clean Benches



WORLD CLASS. WORLDWIDE.

Esco Technologies, Inc. • 2940 Turnpike Drive, Units 15-16 • Hatboro, PA 19040, USA  
Toll-Free USA and Canada 877-375-ESCO • Tel 215-441-9661 • Fax 215-441-9660  
[us.escoglobal.com](http://us.escoglobal.com) • [usa@escoglobal.com](mailto:usa@escoglobal.com)

Esco Micro Pte. Ltd. • 21 Changi South Street 1 • Singapore 486 777  
Tel +65 6542 0833 • Fax +65 6542 6920 • [mail@escoglobal.com](mailto:mail@escoglobal.com)  
[www.escoglobal.com](http://www.escoglobal.com)

Esco Global Offices | Singapore | Philadelphia, USA | Leiden, The Netherlands | Salisbury, UK  
Kuala Lumpur, Malaysia | Beijing, Shanghai & Guangzhou, China | Mumbai, India | Manama, Bahrain

